


## STREPTAVIDIN COATED 12x75 mm TUBES

### Binding capacity streptavidin coated 12x75 mm tubes

Streptavidin coated tubes 12x75 mm were incubated with biotin solutions (from 0 to 195.2 ng/ml) containing 4 ng/ml of biotinylated peroxidase for 30' R.T. After washing step, the streptavidin tubes 12x75 were incubated with TMB and blocked with sulphuric acid 1N. The O.D. values were read at 450 nm.

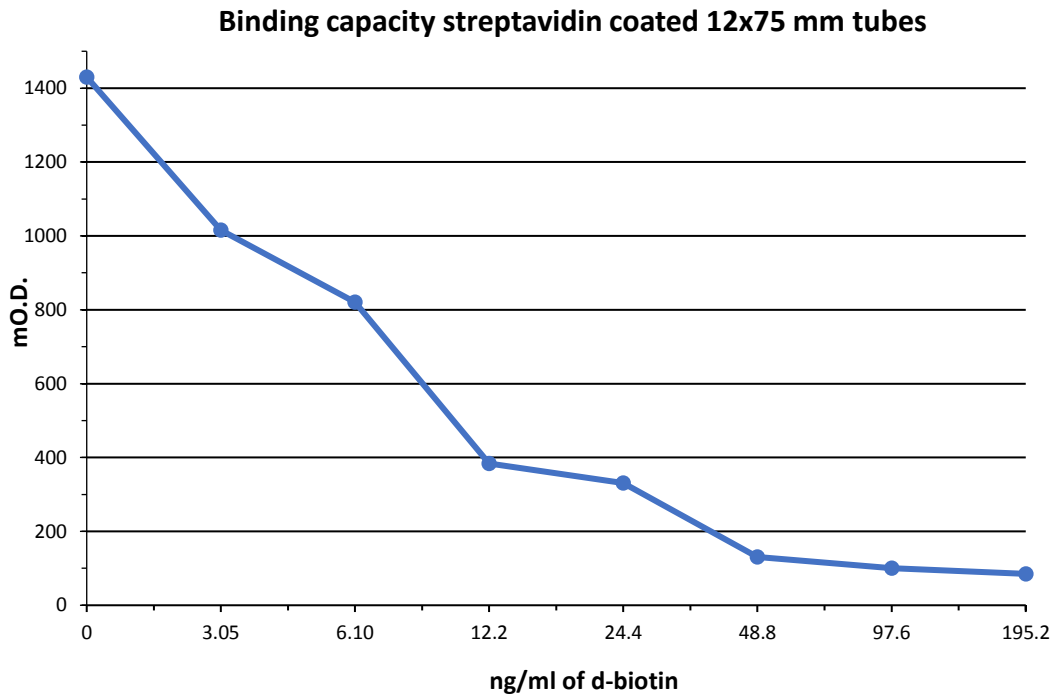

The Biomat Streptavidin 12x75 tubes show a nominal **binding capacity of ~ 28 - 30 pmol d-biotin/tube**

uniformity	CV% < 5
------------	---------