


## STREPTAVIDIN COATED PCR 8 STRIP TUBES

### Binding capacity streptavidin coated PCR strip tubes

Streptavidin coated PCR 8 strip tubes were incubated with biotin solutions (from 0 to 195.2 ng/ml) containing 1.3 ng/ml of biotinylated peroxidase for 30' R.T. After washing step, the PCR tubes strips were incubated with TMB and blocked with sulphuric acid 1N. The O.D. values were read at 450 nm.

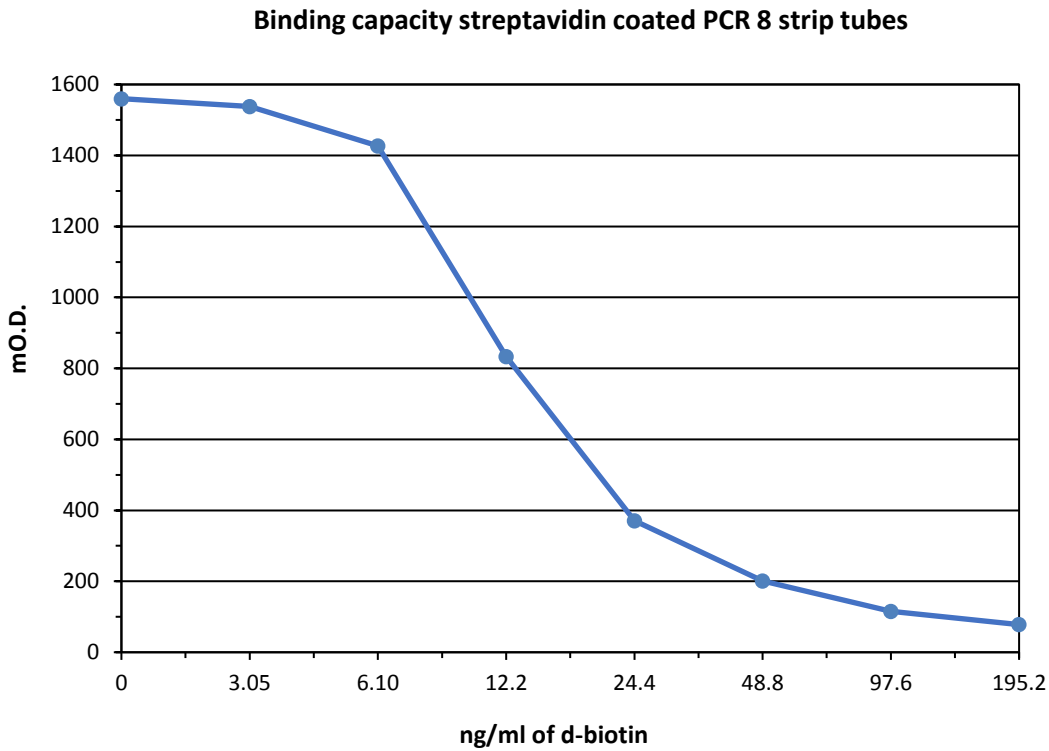

The Biomat Streptavidin PCR 8 tubes strips show a nominal **binding capacity of ~ 5 - 6 pmol d-biotin/tube**