


STREPTAVIDIN HB COATED 12x75 mm TUBES

Binding capacity streptavidin HB coated 12x75 tubes

Streptavidin HB coated tubes 12x75 mm were incubated with biotin solutions (from 0 to 195.2 ng/ml) containing 4 ng/ml of biotinylated peroxidase for 30' R.T. After the washing step, the streptavidin HB tubes 12x75 were incubated with TMB and blocked with sulphuric acid 1N. The O.D. values were read at 450 nm.

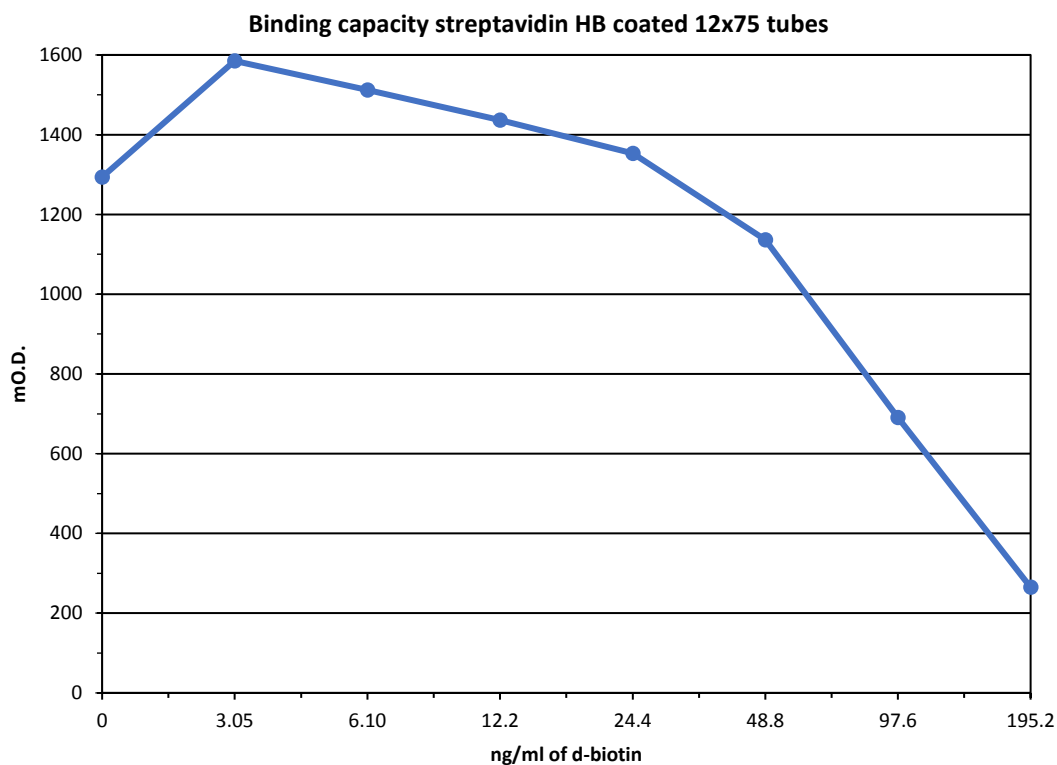

The Biomat Streptavidin 12x75 mm tubes show a nominal **binding capacity of > 400 pmol d-biotin/tube**

uniformity	CV% < 5
------------	---------